

Alternative Methods to RSK 175 Using Purge and Trap Concentration and Automated Headspace for the Analysis of Dissolved Gases in Drinking Water

February 25, 2013

Presenter:

Nathan Valentine, Applications Manager



TELEDYNE TEKMAR
Everywhereyoulook™

Due to increased concern over the hydraulic fracturing process and the release of methane and other chemicals into the local drinking water, a need has developed for fast and accurate analysis of dissolved gasses in water.



TELEDYNE TEKMAR
Everywhereyoulook™

SOP RSK-175

“Sample Preparation and Calculations for Dissolved Gas Analysis in Water Samples Using a GC Headspace Equilibration Technique”

Not an official EPA-approved Method

Analytes: hydrogen, methane, ethylene, ethane, propane, butane, acetylene, nitrogen, nitrous oxide and oxygen

No standard calibration prep method – varies lab to lab

RSK 175 has been employed for analysis of light hydrocarbons in drinking water surrounding hydraulic fracturing well sites



TELEDYNE TEKMAR
Everywhereyoulook™

BOL 6019 (PA-DEP 3686)

Modified method, developed by the Pennsylvania Department of Environmental Protection (PADEP), and helps to simplify and standardize the sample preparation associated with dissolved gas analysis

- Automated headspace analysis
- Liquid calibrations using saturated solutions
- Requires sample handling – VOA => HS vial

Calibration Requirements: %RSD \leq 20% or $r^2 \geq 0.995$

Still requires sample manipulation for each analysis

Limited by equilibration time and platen positions of headspace analyzer



RESULTS

| Compound | Calibration Range | Linearity (r^2) | | MDL | |
|----------|--------------------|---------------------|-------|--------|--------|
| | | HT3 | Versa | HT3 | Versa |
| Methane | 20 ppb to 24 ppm | 0.998 | 0.996 | 2 ppb | 2 ppb |
| Ethene | 141 ppb to 169 ppm | 0.998 | 0.997 | 25 ppb | 19 ppb |
| Ethane | 66 ppb to 79 ppm | 0.999 | 0.999 | 5 ppb | 7 ppb |
| Propane | 74 ppb to 88 ppm | 0.999 | 0.999 | 6 ppb | 4 ppb |



CONCLUSIONS

Both the HT3 and Versa Headspace Analyzers automate the RSK 175 analysis, allowing for higher efficiency and throughput

- **HT3 multi-position platen allows for continuous sample prep**

Decreases potential variability associated with manual GC injections

Met all method performance criteria for all constituents

Standardization to liquid calibrations provides apples to apples comparison to real world samples



TELEDYNE TEKMAR
Everywhere you look™

CAN WE USE PURGE AND TRAP FOR RSK 175?

Trapping Material – Can we trap light gases?

Dynamic Range – low ppb to ppm levels

Purge and Trap Parameters

- Sample Volume
- Purge Volume – Time and Flow

Autosampler

- Sample Temperature
- Carryover

GC Considerations

- Column
- Oven Program
- Split?



TELEDYNE TEKMAR
Everywhere you look™

RESULTS

| Compound | Calibration Range | Linearity (r^2) | %RSD | MDL | % Carryover |
|----------|----------------------|---------------------|------|---------|-------------|
| Methane | 7.92 ppb to 19.8 ppm | 1.000 | 2.0 | 0.4 ppb | 0.04% |
| Ethene | 56.2 ppb to 281 ppm | 0.9995 | 4.5 | 31 ppb | 0.03% |
| Ethane | 26.4 ppb to 132 ppm | 0.9998 | 13.9 | 21 ppb | 0.04% |
| Propane | 29.4 ppb to 147 ppm | 0.9999 | 12.0 | 18 ppb | 0.04% |



CONCLUSIONS

Utilizes existing P&T instrumentation with limited modifications

- Trap, column, and recirculating bath are all that is required

Same advantages to running automated headspace, but also self-contained in 40 mL VOA vials

- No need to manipulate the samples

Potential for fastest run times and highest throughput of all available RSK 175 testing methods since there is no equilibration time

- GC cycle is the limiting factor: too fast = coelutions

New Method: PA-DEP 9423 (October 2012)



TELEDYNE TEKMAR
Everywhereyoulook™